


We provide information and advice to our customers to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose. ChromoCult® is a registered trademark of Merck KGaA, Darmstadt, Germany.



W 286108 03/09

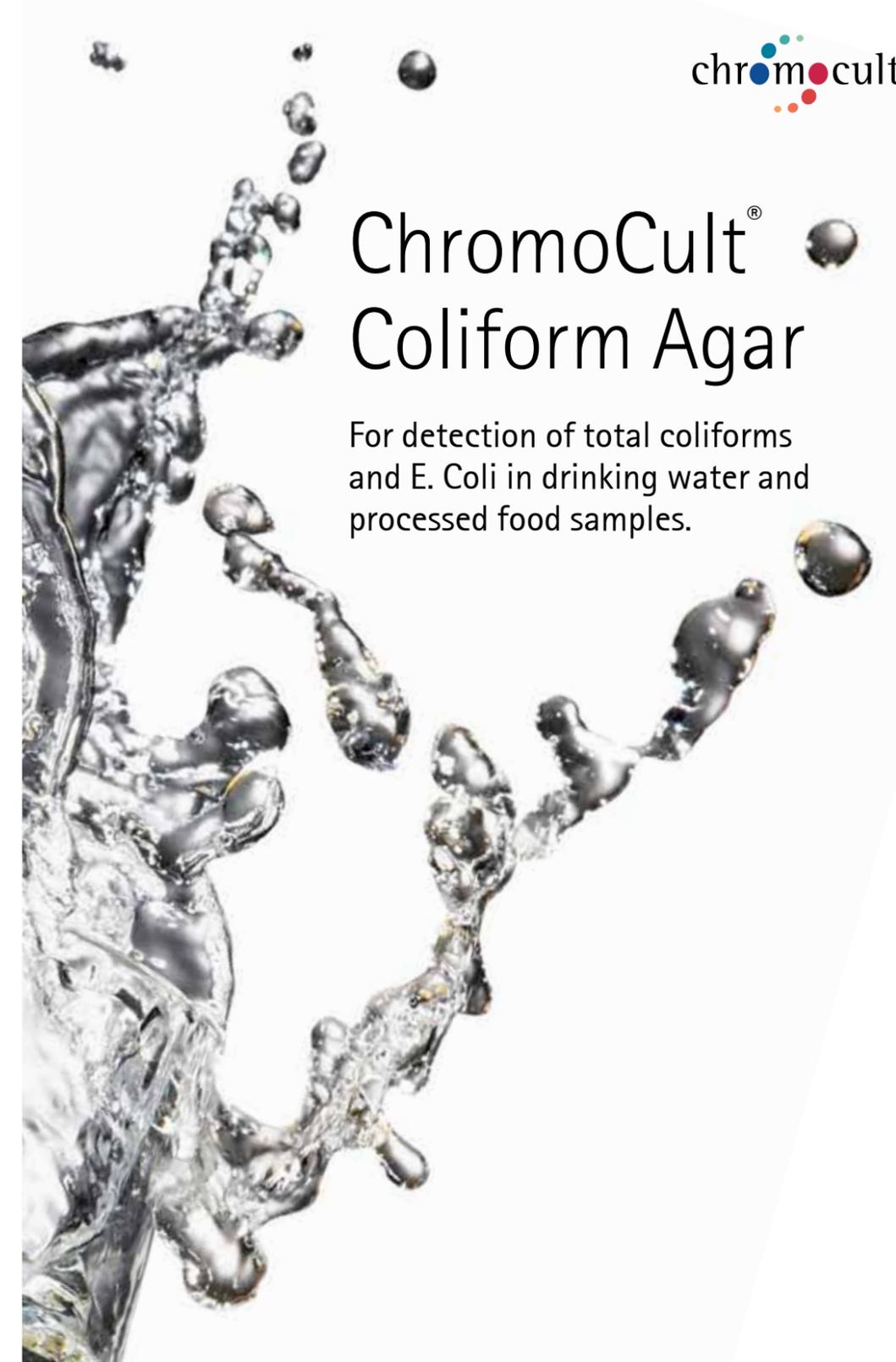

chromocult


# ChromoCult® Coliform Agar

For detection of total coliforms  
and E. Coli in drinking water and  
processed food samples.

Merck KGaA  
64271 Darmstadt, Germany  
Fax +49 (0)61 51/72-60 80  
E-mail: [mibio@merck.de](mailto:mibio@merck.de)  
Internet: [www.merck-chemicals.com](http://www.merck-chemicals.com)



# ChromoCult® Coliform Agar

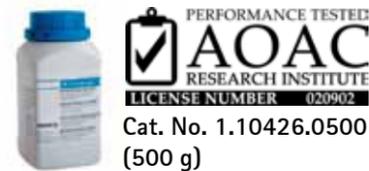

Colour makes the difference.

Selective agar for the simultaneous detection of total coliforms and *E. coli* in drinking water and processed food samples.  
US-EPA approved as alternative method for detection of total coliforms and *E. coli* in Water.  
EU approved as alternative method for detection of *E. coli* and coliforms (in Spain).

The manufacturing process for food and water often result in metabolic injury to the microorganisms and/or reduction of microbial members. Sub lethal microbial injury is caused by chemical or heat treatment as well as storage under frozen, dried or chilled conditions.

Traditional culture media used for the detection of coliforms/*E. coli* often contain ox bile or bile salts to inhibit Gram-positive bacteria. These strong inhibitors however may limit the growth of the targeted microorganisms if they are already sub lethally damaged.

## Intended Use

ChromoCult® Coliform Agar is a selective and differential chromogenic culture medium intended for use in microbiology laboratories analyzing food and water. Within 24 hours this medium enables the detection, differentiation and enumeration of *E. coli* and coliforms from drinking water and processed food matrices such as frankfurters, cooked chicken and non-fat dried milk.

A careful selection of inhibitors used in the selective media is required to ensure the growth and recovery of these damaged microorganisms.

ChromoCult® Coliform Agar contains Tergitol® 7 as an inhibitor of Gram-positive bacteria which has no negative effect on the growth of the targeted coliforms/*E. coli*. ChromoCult® Coliform Agar is therefore the ideal medium for the detection of coliforms/*E. coli* in drinking water and processed foods.

## Application

### Water testing

For water testing the agar medium base should be used with the addition of antibiotics. In the USEPA study the agar medium base was supplemented with 10 mg of cefsulodin per liter. The European study in Spain was performed with the agar medium base supplemented with the *E. coli*/Coliform Selective Supplement. ChromoCult® Coliform Agar is usually inoculated by the membrane filter method.

Filter an appropriate volume of sample (e.g., 100 ml for municipal drinking water, 250 ml for bottled water, or 1 ml for surface water) using a membrane filter. Place the filter on ChromoCult®

Coliform Agar ensuring that no air is trapped underneath. Incubate the inoculated dishes aerobically at 35–37 °C in an inverted position for 24 hours.

After incubation, examine the plates for the presence of typical colonies of *E. coli* and other coliforms.

NOTE: The type of membrane filter affects the size and coloration of colonies. The best results were obtained using membrane filters of cellulose mixed-ester material, e.g. Pall Gelman GN-6 (OSS-MER et al., 1999).

### Food testing

For processed food samples the agar medium base can be used without the addition of antibiotic supplements (as in the AOAC™ study). For fresh food samples with a higher microbial load, ChromoCult® Coliform Agar ES (EMD cat. no. 1.00850) is recommended.

ChromoCult® Coliform Agar is usually inoculated by the pour plate method. Using a sterile pipette, transfer 1 ml of the liquid test sample (or 1 ml from the appropriate dilution) to a sterile Petri dish. Pour into each Petri dish approximately 15 ml of the ChromoCult Coliform Agar while the medium is still liquefied, but only after the medium has been cooled to 45–50 °C in a water bath. Carefully swirl the plate until the inoculum is thoroughly mixed with the medium. Allow the mixture to solidify on a cool horizontal surface.

Incubate the inoculated dishes aerobically at 35–37 °C in an inverted position (agar side up) for 24 hours. After incubation, examine the plates for the presence of typical colonies of *E. coli* and other coliforms.

## Benefits

### Safe time

Fastest method for simultaneous detection of coliforms and *E. coli* in water and processed food. EPA and AOAC approved

### Safe analytic results

Safe detection of injured bacteria because of careful selected inhibitors. Less danger of false negative results

### Easy

Clear detection of coliforms and *E. coli* even in presence high amount of accompanying flora. Just add the recommend Supplement


### Safe money

Easiest confirmation of *E. coli*. Just add a drop of Indole reagent to a colony and the cherry colour shows you whether it is *E. coli* or not.

More detailed Information about validation studies, limitation or performance characteristics are described in the Technical Information.

## Quality control

Test strains	Inoculum (CFU/plate)	% Recovery rate	Colony colour	Salmon GAL	X-Glucuronide
<i>E. coli</i> ATCC 11775	10–100	≥ 70	dark blue to violet	+	+
<i>E. coli</i> DSM 502	10–100	≥ 70	blue to violet	+	+
<i>Citrobacter freundii</i> ATCC 8090	10–100	≥ 70	salmon to red	+	-
<i>Enterobacter aerogenes</i> ATCC 13408	10–100	≥ 70	salmon to red	+	-
<i>Klebsiella pneumoniae</i> ATCC 13883	10–100	≥ 70	salmon to red	+	-
<i>Salmonella enteritidis</i> ATCC 13076	10–100	not limited	colorless	-	-
<i>Enterococcus faecalis</i> ATCC 19433	1000–2000	≤ 0.01			
<i>Bacillus cereus</i> ATCC 11778	1000–2000	≤ 0.01			

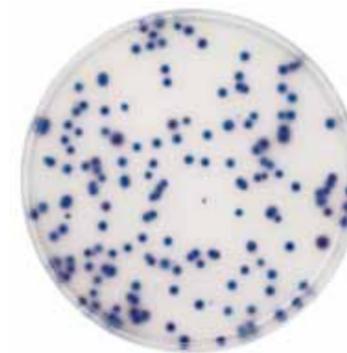

*Escherichia coli* ATCC 11775



*Citrobacter freundii* ATCC 8090

## Ordering Information

Product	Ordering No.	Pack Size
ChromoCult® Coliform Agar	1.10426.0500	500 g
<i>E. coli</i> /Coliform Selective-Supplement	1.00898.0001	1 x 16 vials
Buffered Peptone water (BPW)	1.07228.0500	500 g
Sodium chloride peptone broth (buffered)	1.10582.0500	500 g
Chromoplate® Coliform Agar	1.10156.0001	20 plates

## Literature

- American Public Health Association. 1995. *Standard methods for the examination of water and wastewater*. 19th ed. American Public Health Association, Washington, D.C., USA.
- Byamukama, D., Mach, R.L., Kansime, F., Manafi, M., and A.H. Farnleitner. 2005. *Discrimination Efficacy of Fecal Pollution Detection in Different Aquatic Habitats of a High-Altitude Tropical Country, Using Presumptive Coliforms, Escherichia coli, and Clostridium perfringens Spores*. *Appl. Environ. Microbiol.* 71: 65–71.
- Geissler, K., Manafi, M., Amoros, I., and J.L. Alonso. 2000. *Quantitative determination of total coliforms and Escherichia coli in marine waters with chromogenic and fluorogenic media*. *J. Appl. Microbiol.* 88: 280–285.
- Gonzales, R.D., Tamagnini, L.M., Olmos, P.D., and G.B. de Sousa. 2003. *Evaluation of a chromogenic medium for total coliforms and Escherichia coli determination in ready-to-eat foods*. *Food Microbiology* 20: 601–604.
- ISO International Standardisation Organisation. *Microbiology of food and animal feeding stuffs – General rules for microbiological examinations*. ISO 7218:1996/Amendment 1:2001
- ISO International Standardisation Organisation. *Microbiology of food and animal feeding stuffs – Preparation of test samples, initial suspension and decimal dilutions for microbiological examination. Part 1: General rules for the preparation of the initial suspension and decimal dilutions*. ISO 6887–1:1999.
- ISO International Standardisation Organisation. *Water quality – Sampling. Part 1: Guidance on the design of sampling programmes and sampling techniques*. ISO 5667–1:2006.
- ISO International Standardisation Organisation. *Water quality – General guidance on the enumeration of micro-organisms by culture*. ISO 8199:2005.
- KILIAN, M. and BÜLOW, P. *Rapid diagnosis of Enterobacteriaceae. I. Detection of bacterial glycosidases*. *Acta Pathol. Microbiol. Scand. Sect. B* 84: 245–251 (1976)
- New Zealand Dairy Industry. *Microbiological Methods Manual, Section 48: Product Test Methods – Enteric Indicator Organisms*. NZTM 2; 48.5.1–48.5.10 (1998)
- Ogden, I.D., Brown, G.C., Gallacher, S., Garthwaite, P.H., Gennari, M., Gonzalez, M.P., Jorgensen, L.B., Lunestad, B.T., MacRae, M., Nunes, M.C., Petersen, A.C., Rosnes, J.T., and J. Vliegthart. 1998. *An interlaboratory study to find an alternative to the MPN technique for enumerating Escherichia coli in shellfish*. *International Journal of Food Microbiology* 40: 57–64
- Ossmer, R., Schmidt, W., Mende, U. 1999. *ChromoCult® Coliform Agar – Influence of Membrane Filter Quality on Performance*. – XVII Congreso de la Sociedad, Granada, Spain.
- Suwansonthichai, S., and S. Rengpipat. 2003. *Enumeration of coliforms and Escherichia coli in frozen black tiger shrimp Penaeus monodon by conventional and rapid methods*. *International Journal of Food Microbiology* 81: 113–121.
- Turner, K.M., Restaino, L., and Frampton, E.W. 2000. *Efficacy of ChromoCult Coliform Agar for Coliform and Escherichia coli Detection in Foods*. *J. of Food Prot.* 63: 539–541.
- US-EPA approved method. *Federal Register/Vol. 67, No. 209, 29. October 2002, Rules and Regulations*
- U.S. Food and Drug Administration. 2003. *Bacteriological Analytical Manual Online, Chapter 1: Food Sampling and Preparation of Sample Homogenate*, AOAC International, Gaithersburg, MD., USA.

## Additives

Cat. No.	Product	Pack size
1.09293.0100	KOVACS® Indole reagent	100 ml
1.11350.0001	Bactident® Indole	30 ml
Pall 66278	Cellulose-mixed-ester-GN6 membrane filters	
Schleicher & Schüll 406870	Cellulose-mixed-ester-ME 25/21	

